



## Time course of a trial

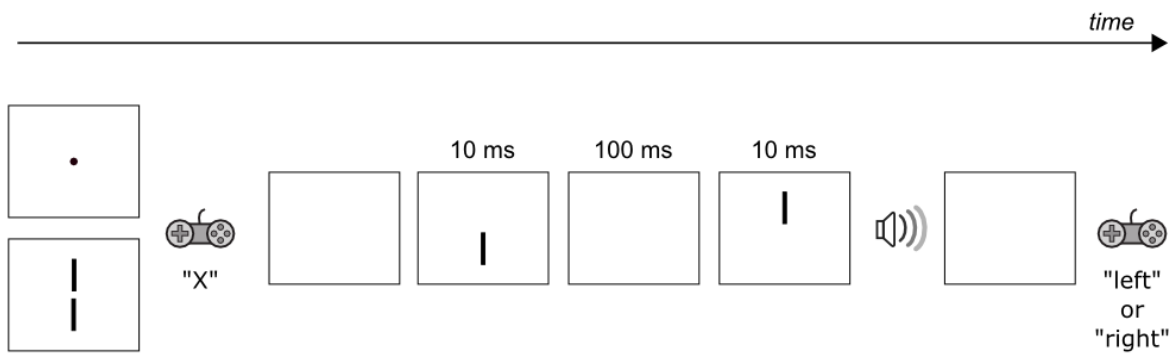


## Task

**Is the top line to the left or to the right of the bottom line?**

## Response

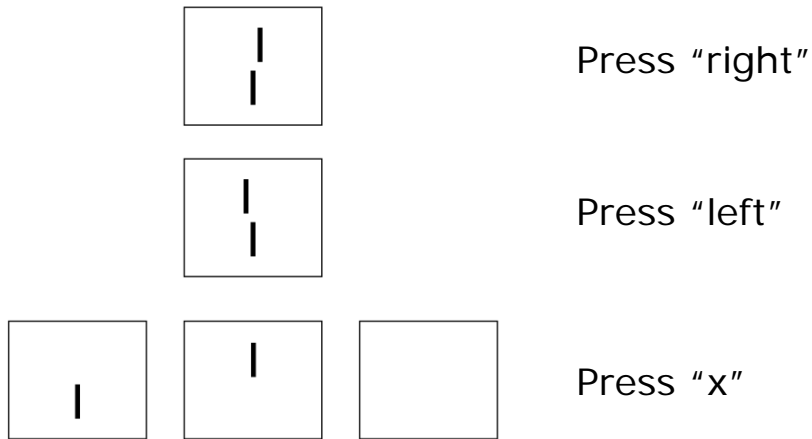


## Stimulus deflector

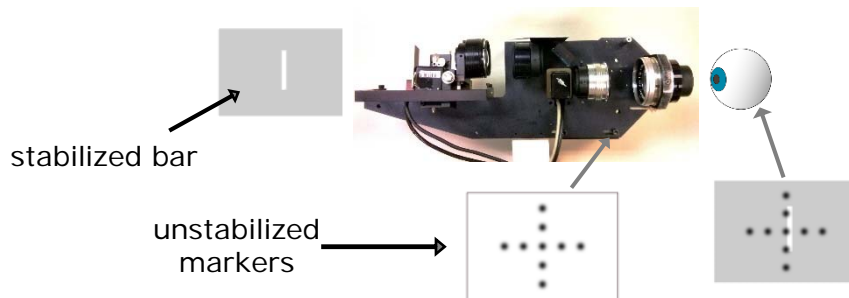

# Calibration





gaze

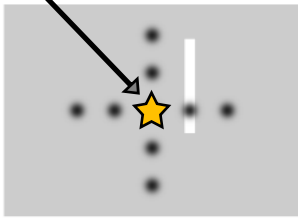



fixate the central unstabilized marker

adjust the **OFFSET** to move the bar at the center of the marker

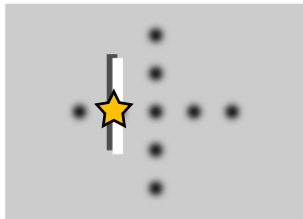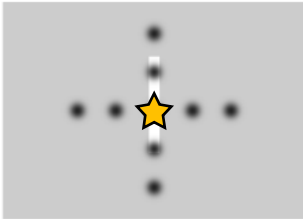


fixate a marker on the periphery on the left

adjust the **AMPLITUDE** to move the bar at the center of the marker (the black bar will disappear)

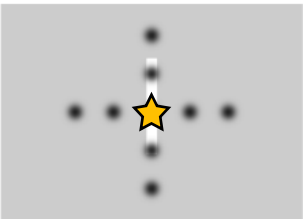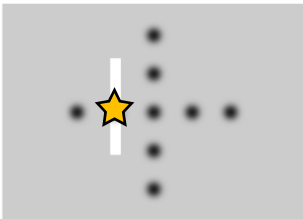


fixate the central unstabilized marker

adjust the **OFFSET** to move the bar at the center of the marker

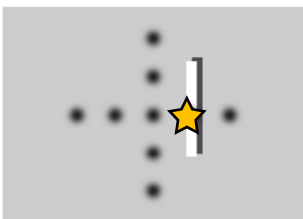

fixate a marker on the periphery on the right

adjust the **AMPLITUDE** to move the bar at the center of the marker (the black bar will disappear)

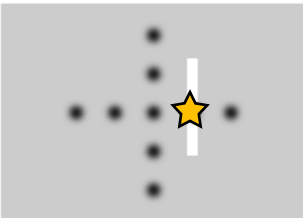

fixate the central unstabilized marker

adjust the **OFFSET** to move the bar at the center of the marker